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NUMERO DE SERIE: _____

- 7 -

HIGH-PRECISION
ELECTRONIC SCALE
LAB-500

USER'S MANUAL
RELEASE: 1

IMPORTANT:
BEFORE OPERATING YOUR SCALE, PLEASE READ THIS MANUAL CAREFULLY.

COD. 13100718


BOX CONTENTS

WITH YOUR PURCHASE YOU WILL RECEIVE:

INSTALLATION INSTRUCTIONS

1.- PLACE THE BATTERIES IN THE BATTERY COMPARTMENT LOCATED UNDER THE SCALE. SLIDE UP THE COVER AND REMOVE IT (USE THE TWO AA BATTERIES INCLUDED)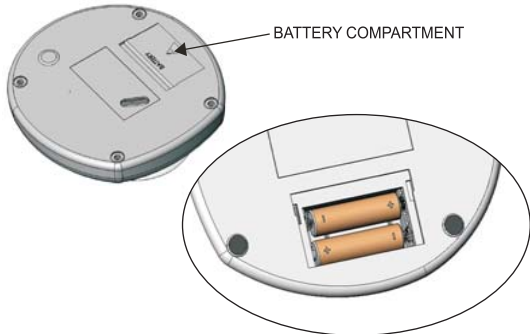

- 1 -

2.- SET THE PLATE OVER THE SCALE'S PLATE-HOLDER BY TURNING IT SMOOTHLY, IN CLOCKWISE DIRECTION, UNTIL IT STOPS.


3.- MAKE SURE THE SCALE IS ON A LEVELLED, VIBRATION-FREE SURFACE.

4.- TURN ON THE SCALE AND WAIT UNTIL THE DISPLAY SHOWS ZERO WEIGHT, THAT MEANS IT IS READY TO BE USED.

OPERATION

THE LAB-500 ELECTRONIC SCALE IS AN EASY-TO-USE PRODUCT. IT CAN BE USED IN LABORATORIES, JEWELRIES, PLASTIC INJECTION, AS WELL AS IN PREPARATION OF PAINT MIX AND MANY OTHER WEIGHING APPLICATIONS. SWITCH BETWEEN GRAMS (gram), POUNDS (Lb), OUNCE (oz) OR CARATS (carat). USE THE TARE FUNCTION TO AUTOMATICALLY SUBSTRACT THE WEIGHT OF A CONTAINER.

KNOWING THE KEYBOARD

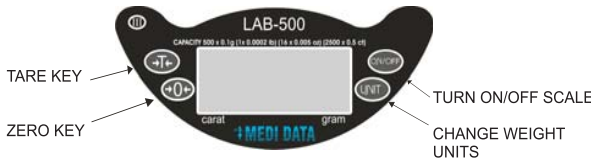- 2 -

HOW TO WEIGH A PRODUCT

PUT THE SCALE IN A FLAT SURFACE AND TURN IT ON BY PRESSING THE **ON/OFF** KEY.

WAIT UNTIL THE ZERO CONDITION IS SHOWN, THEN PLACE THE LOAD ON THE SCALE. SEE THE WEIGHT VALUE ON THE DISPLAY WITH THE DESIRED UNIT BY PRESSING THE "UNIT" KEY. SWITCH BETWEEN "GRAM", "CARAT", "LB" AND "OZ". FOR EXAMPLE, IF A WEIGHT OF 0.5 LB IS PLACED ON THE SCALE, THE DISPLAY WILL SHOW:

HOW TO USE "TARE" FUNCTION

WHEN YOU WANT TO WEIGH A PRODUCT IN A CONTAINER AND DO NOT WANT TO CONSIDER THE WEIGHT OF SAID CONTAINER DURING AN OPERATION, YOU CAN USE THE TARE FUNCTION. TO USE THIS FUNCTION FOLLOW THE FOLLOWING STEPS: SUPPOSE YOU WANT TO WEIGH 0.1 POUNDS OF GOLD DUST BOWL USING THE TARE-PLATE.

1.- PLACE THE EMPTY TARE-PLATE ON THE SCALE.

2.- THE SCALE WILL SHOW THE WEIGHT OF THE TARE-PLATE.

3.- PUSH THE **←T→** KEY. THE DISPLAY WILL SHOW ZERO AND WILL ACTIVE THE **NET** SYMBOL.

4.- PLACE THE PRODUCT IN THE TARE-PLATE. THE SCALE WILL SHOW ONLY THE WEIGHT OF THE PRODUCT WITHOUT THE WEIGHT OF THE TARE-PLATE. NOTE: WHEN THE TARE-PLATE IS REMOVED A NEGATIVE NUMBER (THE WEIGHT OF TARE-PLATE) WILL BE SHOWN IN THE DISPLAY.

- 3 -

IT IS IMPORTANT TO MENTION THAT THIS FUNCTION IS SUBSTRUCTIVE. FOR EXAMPLE IF YOU HAVE A CONTAINER THAT WEIGHS 0.4 POUNDS, THE SCALE WILL ONLY BE ABLE TO WEIGHT UP TO 0.6 POUNDS OF PRODUCT AND COMPLETE THE MAXIMUM CAPACITY (1 POUNDS, 500 GRAMS, 16 OZ OR 2500 CARATS). TO CANCEL THE TARE FUNCTION REMOVE ALL THE WEIGHT OVER THE PLATE AND PRESS THE **←T→** KEY, THE SCALE WILL RETURN TO ITS NORMAL FUNCTION. THE NET SYMBOL WILL DISAPPEAR.

READ THE FOLLOWING OBSERVATIONS CAREFULLY:

A) WHEN THERE IS NO WEIGHT ON THE SCALE YOU CAN NOT ACTIVATE THE "TARE" FUNCTION.

B) WHEN USING THE TARE FUNCTION AND REMOVING THE CONTAINER YOU WILL SEE THE NEGATIVE WEIGHT OF THE CONTAINER. IF YOU REMOVE A LOAD GREATER THAN THE CONTAINER WEIGHT YOU WILL GET A "NEG" MESSAGE

IF IT IS THE CASE, JUST TURN OFF THE SCALE AND TURN IT BACK ON TO GET A NEW ZERO REFERENCE.

HOW TO USE THE **UNIT** KEY

IF THE SCALE IS IN "gram" MODE AND YOU PLACE A PRODUCT WITH A WEIGHT OF 300 GRAMS ON THE PLATE, YOU WILL SEE THE FOLLOWING DIGITS IN THE DISPLAY:

IF YOU PRESS THE **UNIT** KEY, THE SCALE WILL CHANGE TO "carat" MODE AND YOU WILL SEE THE FOLLOWING DIGITS IN THE DISPLAY:

IF YOU PRESS THE **UNIT** KEY AGAIN, THE SCALE WILL CHANGE TO "Lb" MODE AND, YOU WILL SEE THE FOLLOWING DIGITS IN THE DISPLAY:

- 4 -

HOW TO USE THE "ZERO" REFERENCE

IT IS POSSIBLE THAT THE SCALE SHOWS A NUMBER IN THE DISPLAY WITHOUT HAVING ANY PRODUCT ON THE PLATE DUE TO ROUGH HANDLING OR SOME OTHER FACTORS WHEN THIS HAPPENS PRESS THE **←0→** KEY AND THE DISPLAY WILL RETURN TO "ZERO". THE ZERO SYMBOL WILL APPEAR IN THE DISPLAY.

AUTO POWER-OFF TIMER FUNCTION

THIS FUNCTION AUTOMATICALLY TURNS OFF THE SCALE IF THE SCALE HAS NOT BEEN USED FOR A PERIOD OF 5 MINUTES TO SAVE ENERGY CONSUMPTION

LOW BATTERY INDICATOR

WHEN THE SCALE RUNS OUT OF BATTERY, YOU WILL SEE THE LOW BATTERY SYMBOL ON THE DISPLAY, BUT YOU WILL STILL BE ABLE TO OPERATE THE SCALE FOR A SHORT PERIOD OF TIME.

WHEN THE VOLTAGE OF THE BATTERY IS VERY LOW YOU WILL SEE THE "LOW" MESSAGE AS FOLLOWS:

AND YOU WON'T BE ABLE TO OPERATE THE SCALE UNTIL THE BATTERY HAS BEEN REPLACED.

HELP CONDITION

THIS MESSAGE APPEARS IN THREE DIFFERENT SITUATIONS.

A) WHEN THE SCALE WAS TURNED ON WITH AN EXCESSIVE WEIGHT ON THE PLATE (GREATER OR EQUAL TO 50% OF ITS CAPACITY). TO FIX THIS SITUATION REMOVE THE LOAD FROM THE SCALE AND TURN ON IT AGAIN.

B) WHEN THE SCALE WAS HANDLED ROUGHLY AND/OR GOT HIT DURING ITS USE, CAUSING IRREVERSIBLE DAMAGE ON THE WEIGHT SENSOR.

C) WHEN THE SCALE HAS BEEN LOADED WITH A LOAD GREATER THAN THE MAXIMUM CAPACITY (500 GRAMS)

- 5 -

MAINTENANCE AND CLEANING

ONE OF THE GREAT ADVANTAGES OF BUYING THIS TYPE OF SCALE IS THAT IT REQUIRES MINIMAL MAINTENANCE AND IS VERY SIMPLE TO KEEP IT IN A FAVORABLE STATE.

• **DO NOT OVERLOAD THE SCALE WITH A LOAD GREATER THAN 500 GRAMS, SINCE YOU MAY DAMAGE THE HIGH-PRECISION WEIGHING SENSOR**

• **DO NOT USE WATER BLAST OR CORROSIVE LIQUIDS TO CLEAN THE SCALE. USE A WET CLOTH AND WIPE EACH OF THE COMPONENTS.**

• **WHEN CLEANING THE SCALE MAKE SURE THAT THE SPACE UNDER THE PLATE IS COMPLETELY CLEAN, SO NOTHING OBSTRUCTS THE SCALE OPERATION.**

• **DO NOT EXPOSE THE SCALE TO TEMPERATURES BELOW 32 °F (0 °C) OR ABOVE 104 °F (40 °C).**

• **NEVER PRESS THE KEYBOARD WITH SHARP OBJECTS SUCH AS PENCILS, PENS, RAZORS, ETC.**

• **DO NOT DROP THE PRODUCT TO BE WEIGHED ON THE SCALE PLATE BECAUSE IT WILL DAMAGE THE LOAD CELL.**

• **DO NOT LIFT YOUR SCALE FROM THE PLATE. ALWAYS LIFT IT FROM THE BASE.**

TROUBLESHOOTING GUIDE

PROBLEM	POSSIBLE CAUSE	SOLUTION
IT DOES NOT OPERATE	THE BATTERIES ARE NOT INSTALLED	CHECK THE BATTERY COMPARTMENT
	THE BATTERIES ARE INCORRECTLY INSTALLED	CHECK THAT THE BATTERY TERMINALS(+,-) MATCH WITH THE INDICATION IN THE COMPARTMENT
	LOW BATTERY	REPLACE THE BATTERY
THE SCALE SHOWS DIFFERENT WEIGHTS WITH THE SAME LOAD SEVERAL TIMES	THERE IS MATERIAL UNDER THE PLATE THAT IS OBSTRUCTING THE CORRECT OPERATION OF THE SCALE	REMOVE ANY MATERIAL UNDER THE PLATE. MAKE SURE IT IS VERY CLEAN.


- 6 -

TECHNICAL SPECIFICATIONS

MODEL	LAB-500
MAX. CAPACITY	500 gram, 1 lb, 16 oz, 2500 carat
MIN. DIVISION	0.1 gram, 0.0002 lb, 0.005 oz, 0.5 carat
MAX. TARE	199.9 gram, 0.4408 lb, 7.050 oz, 999.5 carat

OPERATING TEMPERATURE RANGE	32 TO 104 °F (0 TO 40 °C)
STORAGE TEMPERATURE RANGE	14 TO 122 °F (-10 TO 50 °C)

DIMENSIONS (inches)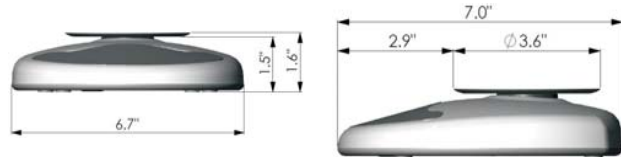

NOTE: THIS PRODUCT SERIAL NUMBER IS LOCATED UNDER THE SCALE. WE SUGGEST YOU TO WRITE DOWN THE SERIAL NUMBER AND MODEL ON THE SPACE BELOW AND KEEP THEM AS PERMANENT PROOF OF YOUR PURCHASE TO HELP IDENTIFY IT IN CASE OF LOSS OR FURTHER ASSISTANCE.

DISTRIBUTOR: _____

ADDRESS: _____

TELEPHONE: _____

DATE OF PURCHASE: _____

MODEL: _____

SERIAL NUMBER: _____

- 7 -